

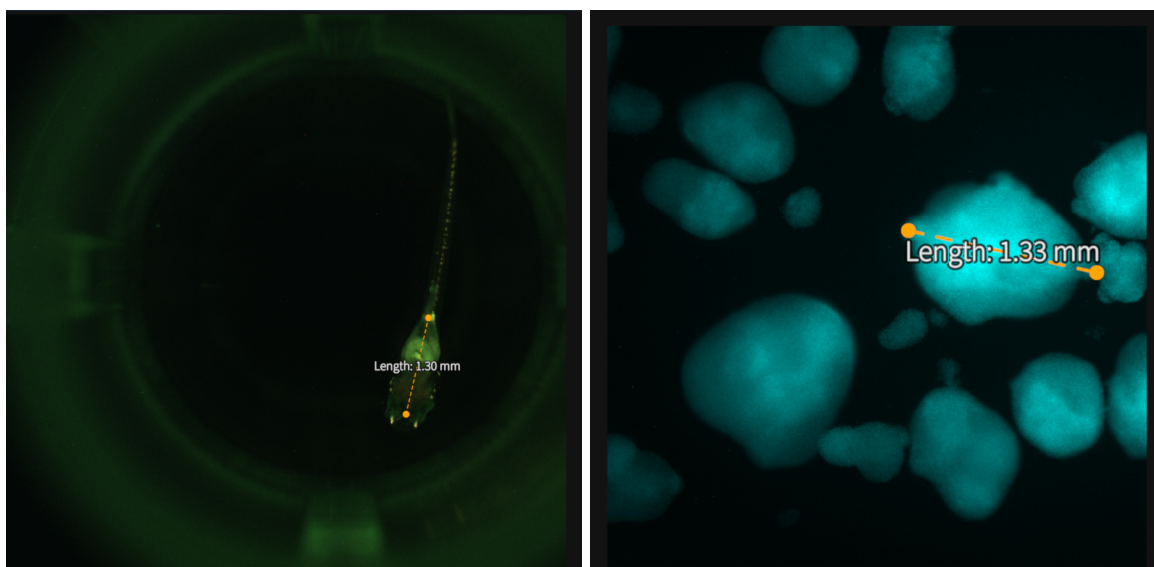


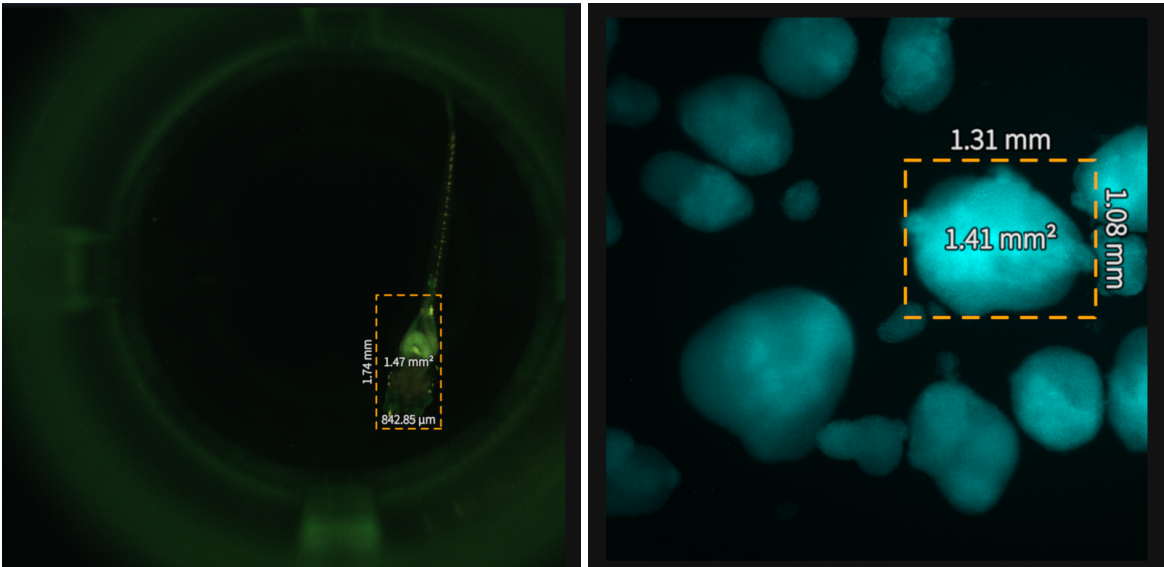

Step	Function
1	Open the MCAM software.
2	Go to Tools > Length Tool.
3	Click once with the left mouse button to define the beginning of the line, move the cursor to extend the line and finally click again with the left mouse button to define the end of the line.
4	To take another measurement simply repeat step 3 . To remove the Length Tool repeat step 2 .

Rectangle Tool

When the Rectangle Tool is selected, rectangles can be drawn across a desired area of the screen. Height, width, area and their corresponding units will be displayed on the screen.

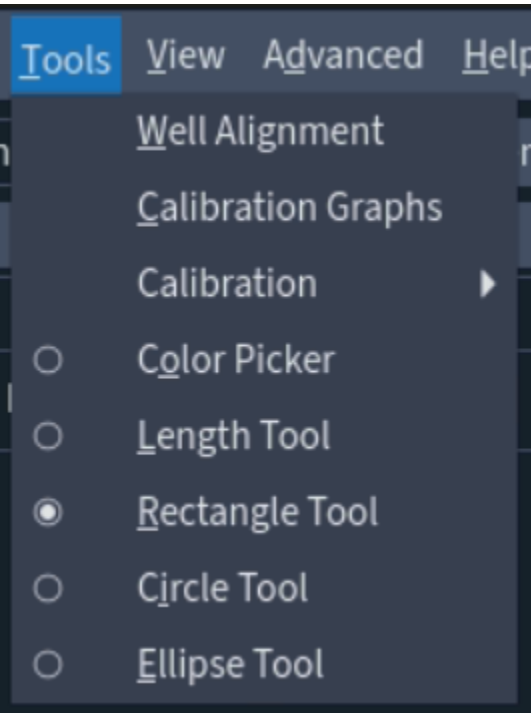

Step	Function
1	Open the MCAM software.
2	Go to Tools > Rectangle Tool.

3	Click once with the left mouse button to define one corner of the rectangle, move the cursor to extend the area covered by the rectangle and click again with the left mouse button to define the second defining corner of the rectangle.
4	To take another measurement simply repeat step 3 . To remove the Rectangle Tool repeat step 2 .

Circle Tool

When the Circle Tool is selected, circles can be drawn across a desired area of the screen. Diameter, area and their corresponding units will be displayed on the screen.

