

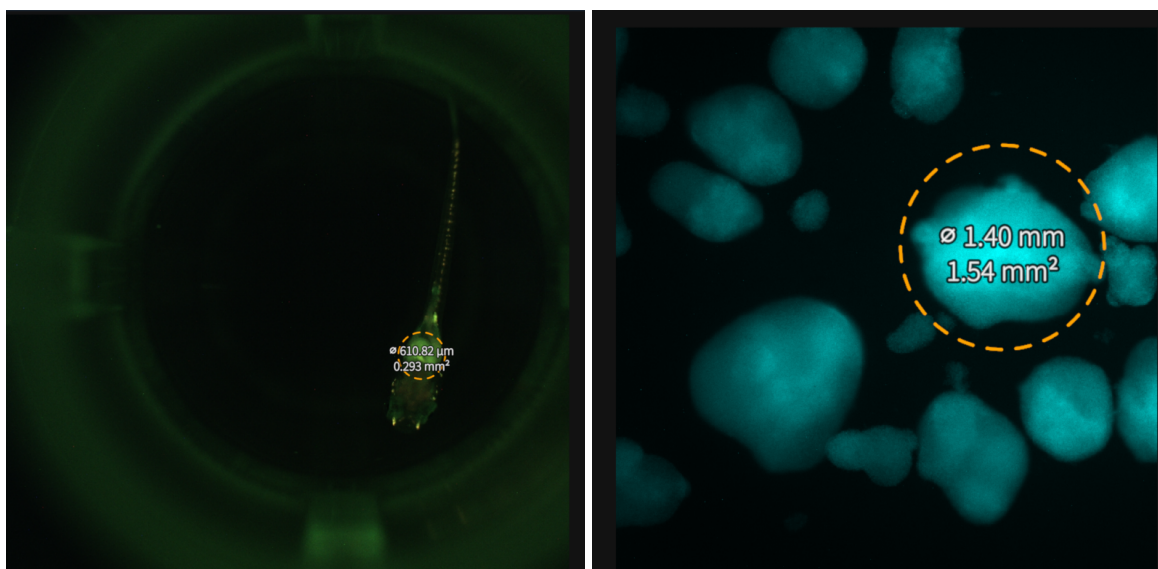


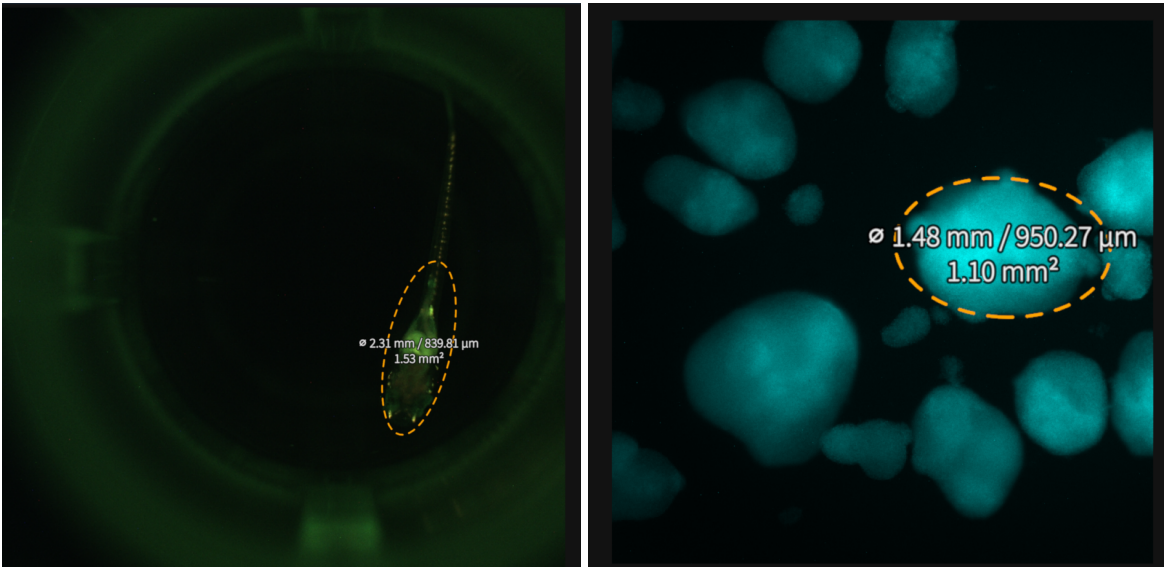

| Step | Function                                                                                                                                                                                                                                       |
|------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| 1    | <b>Open the MCAM software.</b>                                                                                                                                                                                                                 |
| 2    | <b>Go to Tools &gt; Circle Tool.</b>                                                                                                                                                                                                           |
| 3    | <b>Click once with the left mouse button</b> to define the center of the circle, <b>move the cursor</b> to extend the circle's radius and finally click once with the left mouse button a second time to define the final shape of the circle. |
| 4    | To take another measurement simply <b>repeat step 3</b> . To remove the Circle Tool <b>repeat step 2</b> .                                                                                                                                     |

## Ellipse Tool

When the Ellipse Tool is selected, ellipses can be drawn across a desired area of the screen. Width, height, area and their corresponding units will be displayed on the screen.

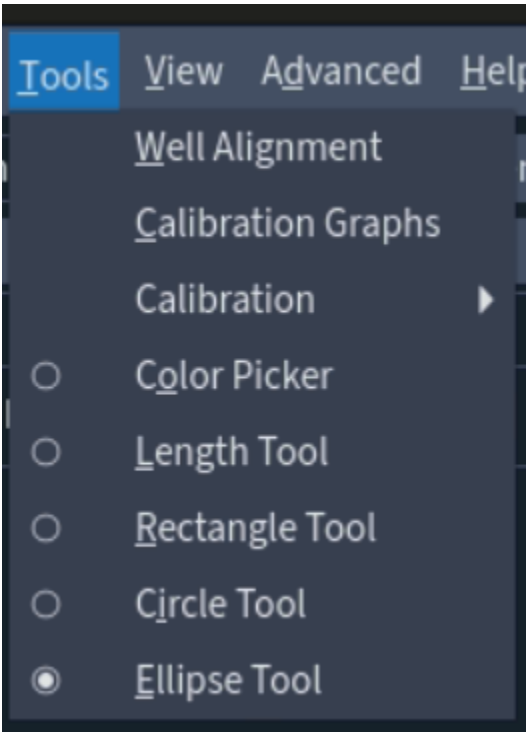

| Step | Function                    |
|------|-----------------------------|
| 1    | Open the MCAM software.     |
| 2    | Go to Tools > Ellipse Tool. |

|   |                                                                                                                                                                                                                                                                                        |
|---|----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| 3 | <b>Start by creating a circle</b> by clicking once with the left mouse button to define the center of the circle. <b>Move the cursor</b> to extend the circle's radius and finally <b>click once with the left mouse button</b> a second time to define the final shape of the circle. |
| 4 | Once the desired circle is created, <b>click once with the left mouse button</b> to shape the ellipse based on this circle. <b>Move the cursor</b> to adjust the ellipse shape/direction and <b>click once with the left mouse button</b> to define the final ellipse shape.           |
| 5 | To take another measurement simply <b>repeat steps 3 and 4</b> . To remove the Ellipse Tool <b>repeat step 2</b> .                                                                                                                                                                     |