



Step	Function
1	Open the MCAM software.
2	Go to Tools > Color Picker.
3	Click on a pixel in the image to get the displayed color at the location.
4	In the MCAM software, go to View which will give the user several options to alter the viewing color space. Changing this will alter the pixel values displayed by the color picker.
5	To take another measurement simply repeat step 3 . To remove the Color Picker repeat step 2 .

Length Tool

When the Length Tool is selected, straight lines can be measured on the screen. The measurement and unit will be displayed on the screen.

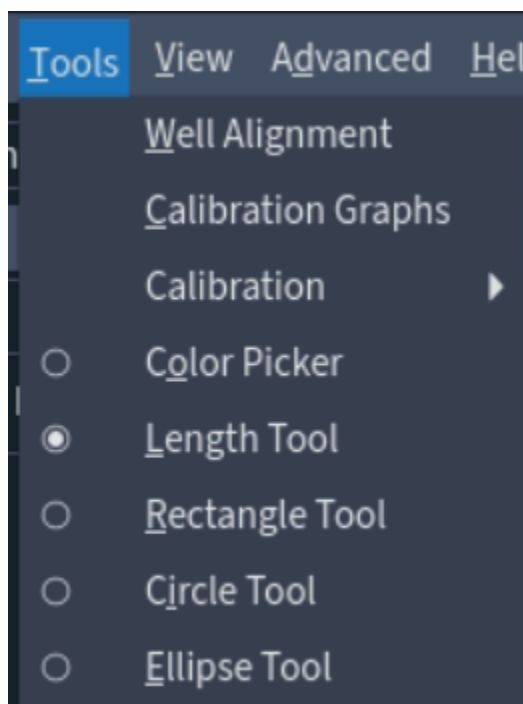

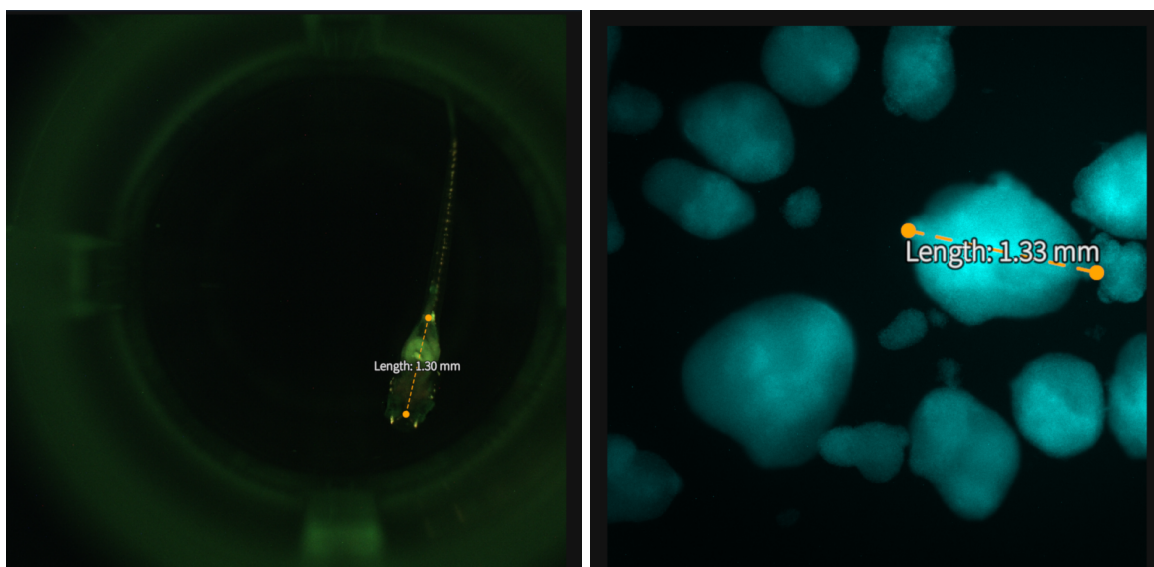

Step	Function
1	Open the MCAM software.
2	Go to Tools > Length Tool.
3	Click once with the left mouse button to define the beginning of the line, move the cursor to extend the line and finally click again with the left mouse button to define the end of the line.
4	To take another measurement simply repeat step 3 . To remove the Length Tool repeat step 2 .

Rectangle Tool

When the Rectangle Tool is selected, rectangles can be drawn across a desired area of the screen. Height, width, area and their corresponding units will be displayed on the screen.