


Acquire an Image

Capturing an image is the simplest form of data acquisition within the MCAM platform.

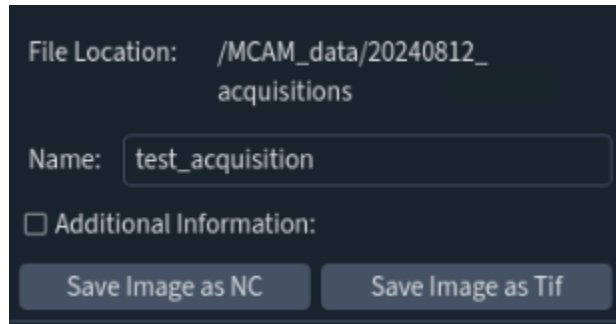

File Location: /MCAM_data/20240812_acquisitions

Name: test_acquisition

☐ Additional Information:

Save Image as NC Save Image as Tif

Step	Function
1	Open the MCAM software.
2	In the left side panel, in the top left corner, click "Select a File Location" . Always select a folder on the "MCAM_data" drive to store data from MCAM acquisitions. The "File Location" will be displayed in the left side panel once defined.
3	Enter a Filename for the image. A timestamp will automatically be appended to the chosen filename.
4	Adjust imaging settings in the left side panel to optimize image quality.
5	Click Save Image as NC or Save Image as Tif to save the image.