



3	Click once with the left mouse button to define one corner of the rectangle, move the cursor to extend the area covered by the rectangle and click again with the left mouse button to define the second defining corner of the rectangle.
4	To take another measurement simply repeat step 3 . To remove the Rectangle Tool repeat step 2 .

Circle Tool

When the Circle Tool is selected, circles can be drawn across a desired area of the screen. Diameter, area and their corresponding units will be displayed on the screen.

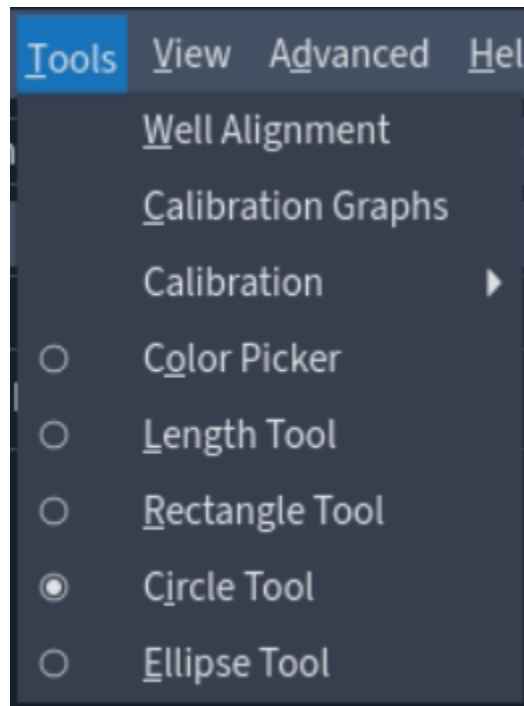

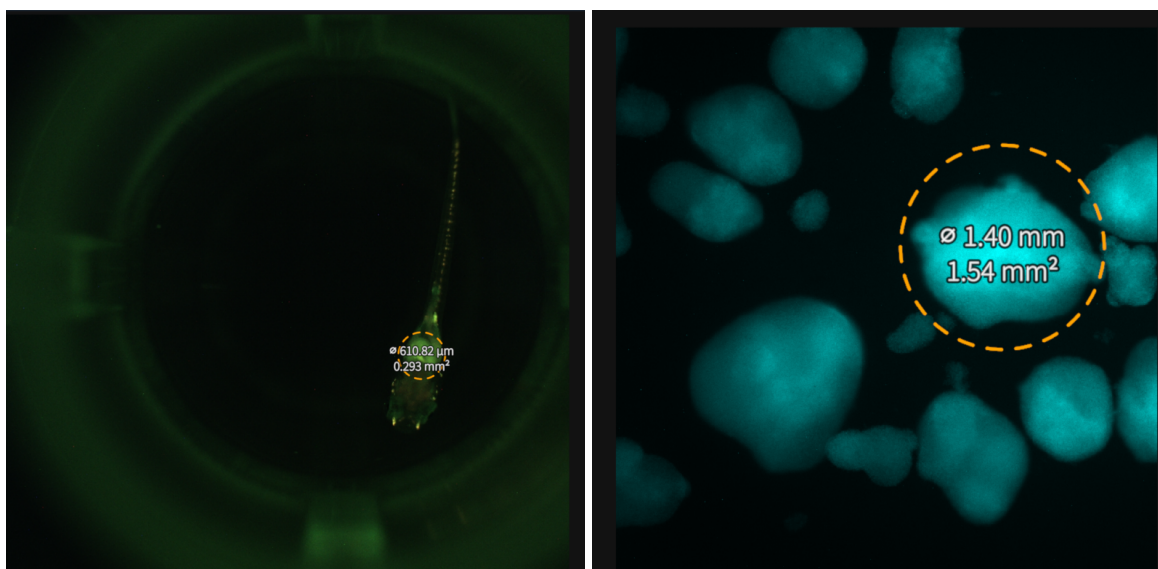

Step	Function
1	Open the MCAM software.
2	Go to Tools > Circle Tool.
3	Click once with the left mouse button to define the center of the circle, move the cursor to extend the circle's radius and finally click once with the left mouse button a second time to define the final shape of the circle.
4	To take another measurement simply repeat step 3 . To remove the Circle Tool repeat step 2 .

Ellipse Tool

When the Ellipse Tool is selected, ellipses can be drawn across a desired area of the screen. Width, height, area and their corresponding units will be displayed on the screen.